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Instruction and maintenance manual
Livret d'instruction et entretien
Bedienungs- und Wartungsanleitung

ARMADIO REFRIGERATO
REFRIGERATED CABINET
ARMOIRE REFRIGEREE
KÜHLSCHRANK

VERS. 1119

ITALIANO

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ENGLISH

Content

Sect. 1 Standards and general warnings	12	4.3 Safety dataplates and guards.....	16
1.1 Testing and guarantee.....	12	4.4 Operating limits	16
1.2 Introduction.....	12	Sect. 5 Routine and programmed maintenance	16
1.3 Product description.....	12	5.1 Basic safety regulations	16
1.4 General safety regulations	13	5.2 Cleaning the refrigerator.....	17
1.5 Client's responsibilities	13	5.3 Periodic checks	17
1.6 Client service requests	13	5.4 Precautionary measures for prolonged disuse.....	17
1.7 Ordering spare parts	13	5.5 Preventive maintenance.....	17
Sect. 2 Specifications	13	Sect. 6 Special maintenance and repairs	18
2.1 Product configuration	13	Sect. 7 Diagnostics	18
2.2 Power output and absorbed power	13	Sect. 9 Symbols	26
2.3 Noise level.....	14		
2.4 Materials and refrigerants.....	14		
Sect. 3 Installation	14		
3.1 Transport and handling	14		
3.2 Positioning	14		
3.3 Wiring and electrical hook-up.....	14		
3.4 Set-up operations.....	15		
3.5 Re-installation.....	15		
3.6 Scrapping and disposal	15		
Sect. 4 Operation	15		
4.1 Applications and intended use	15		
4.2 Safety and accident prevention.....	15		

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Section 1 STANDARDS AND GENERAL WARNINGS

1.1 TESTING AND GUARANTEE

The appliance is tested in our works in compliance with established regulations and then shipped ready for use.

The guarantee is valid for a full 12 months from the date of delivery of the appliance and it covers the repair or replacement of any defective parts, with the exception of electrical and electronic components.

Manifest defects or differences with respect to the client's order must be communicated to the manufacturer within five days from the receipt of the goods or they will not be covered by the guarantee terms.

Any hidden or other defects must be communicated to the manufacturer within five days from the time that they are discovered and, in any event, within the maximum guarantee term of six months. The purchaser shall be entitled only to request repair or replacement of the goods. The purchaser is not entitled to claim compensation for direct or indirect damages of any whatsoever nature. In any event, the entitlement to repair or replacement of the materials must be exercised within the maximum term of the guarantee, which is contractually stipulated to cover a shorter period than the maximum term established by law.

Repairs or replacement of defective materials will be carried out at the manufacturer's works; material returned to the manufacturer must be shipped carriage paid and will be returned to the purchaser carriage forward.

1.2 INTRODUCTION

This manual has been prepared with the scope of supplying all the instructions required for the correct use of the appliance and to maintain it in optimal condition. It also contains important user safety information.

The following professional roles are explained in order to define the responsibilities of each:

Installer: a qualified technician who positions the appliance and places it in service in accordance with the instructions in this manual.

User: the person who, after reading this manual carefully, operates the appliance in accordance with the intended use specified in this manual.

Users' responsibilities:

- to ensure that the products are conserved at suitable temperatures and not exceeding the permitted period of time
- to be aware of the regulations governing the conservation of products and to observe any whatsoever hygiene indications that may be applicable.

The user is obliged to read the manual attentively and refer to the information in the manual at all times.

Particular attention must be paid to the contents of heading 1.5 General Safety Warnings.

Routine Maintenance Technician: qualified technician able to perform routine maintenance of the appliance by following the instructions in this manual (see section 5).

Special Maintenance Technician: qualified technician, authorized by the manufacturer to perform extraordinary maintenance of the appliance (see section 6).

The symbol appears at certain points in the manual to draw the reader's attention to important safety information.

The manufacturer declines any whatsoever responsibility in the case of improper use of the appliance deviating from the reasonably construed intended use, and for all operations carried out that are not in compliance with the instructions laid down in the manual.

This manual must be conserved in a place that is accessible and known to all operators (installer, user, routine maintenance technician, special maintenance technician).

This manual must not be reproduced or divulged, in whole or in part, using any whatsoever means or in any whatsoever form.

1.3 PRODUCT DESCRIPTION

The appliance comprises a modular single body with panelling in various materials and insulation in expanded polyurethane foam, density 42 kg/cu.mt.

The appliance instruments are located on the front panel which in some models close the front of the motor unit, inside which the condenser unit and electrical wiring can be housed.

The appliance interior is fitted with suitable supports for wire shelves (grids) and/or other accessories (drawers and/or baskets)

The doors are fitted with an automatic return device and magnetic seal elements.

During the design and construction stage all measures have been adopted to implement total safety including radiused interior corners, funnel-shaped base panel to convey condensate to exterior, no rough surfaces, fixed guards protecting moving or potentially dangerous parts.

1.4 GENERAL SAFETY REGULATIONS

Read this manual carefully and follow the prescriptions contained herein.

The user assumes full responsibility in the case of operations carried out without observing the instructions in the manual.

Primary general safety regulations:

- do not touch the unit with wet hands and/or feet
- do not use the appliance with bare feet
- do not insert screwdrivers or other pointed objects between guards or moving parts of the appliance
- do not pull the power cord to disconnect the appliance from the electrical mains
- make sure that the appliance is not used by children or unsuitably qualified persons
- before performing any cleaning or maintenance on the appliance disconnect it from the electrical mains by switching of the main switch and extracting the plug
- in the case of faults or malfunctions, switch off the appliance and do not attempt to repair it yourself. All service and repair operations must be performed exclusively by suitably qualified authorized technicians.

1.5 CLIENT'S RESPONSIBILITIES

The customer is required to:

- execute the electrical connection of the appliance
- prepare the place of installation
- provide consumable materials for cleaning
- perform routine maintenance

In the case of power failures or malfunctions do not open the doors and drawers in order to maintain uniform temperature inside the unit. If the problem persists for more than a few hours, move the stored material to a more suitable place.

1.6 CLIENT SERVICE REQUESTS

For all technical problems and any requests for technical service, refer exclusively to your local dealer.

1.7 ORDERING SPARE PARTS

Spare parts orders must be made by consulting the relative spare parts catalogue which gives the correct description of the part, the part reference code and the serial number of your appliance.

The manual is available, upon request, by the manufacturer.

Section 2 SPECIFICATIONS

2.1 PRODUCT CONFIGURATION

The products must be stored in observance of the load limits allowed by the shelves, drawers or baskets in order to ensure efficient air circulation inside the appliance.

DESTINATION OF USE		
MPR	+2°C / +15°C	Biological refrigerators for medical-pharmaceutical use
MPRR	+2°C / +15°C +2°C / +15°C	
LR	0°C / +15°C	Biological refrigerators for laboratory use
LCRR	+0°C / +15°C +0°C / +15°C	
BLFG	-5°C / -25° C	Laboratory freezers
LFG	-5°C / -20° C	
LF	-5°C / -20 ~ -25°C	
LDF	-15°C / -30°C	
PDF	-25°C / -40°C	
LCRF	+0°C / +15°C -5°C / -20 ~ -25°C	Combinations of refrigerators / freezers for laboratory
CI	+2°C / +50°C	Cooled incubators
PI	+22°C	Platelet incubators

2.2 POWER OUTPUT AND ABSORBED POWER

Technical data for power output and absorbed power are shown in the appliance's technical identification label.

Bear in mind the operational limit characteristics as indicated in heading 4.4.

2.3 NOISE LEVEL

The noise level of the appliance is below 70 dB (A), change model-by-model and according to the ambient and usage working conditions.

2.4 MATERIALS AND REFRIGERANTS

Materials in contact or potentially in contact with the stored products are in compliance with the relevant directives. The appliance is designed and built so that the parts in contact with the stored material can be cleaned before each use. The refrigerants utilized comply with established regulations. For more information check the appliance's technical identification label.

Section 3 INSTALLATION

3.1 TRANSPORT AND HANDLING

The appliance must be transported and handled exclusively in a vertical position, in observance of the instructions printed on the packing.

This precaution is necessary to avoid contamination of the refrigerant circuit with compressor lube oil with resulting valve and heat exchanger coil failure and problems starting the electric motor.

The manufacturer accepts no responsibility for problems due to transport executed in conditions other than those specified above.

The accessories supplied with the appliance (runners, wire shelves, basins, drawers, baskets) are supplied in separate packs shipped inside or separately from the unit.

The appliance is secured to a wooden base by means of a screw through each foot and wrapped in polyethylene or packed in a carton, cage or crate.

Refer to heading 3.6 for information on correct disposal of packing material.

The appliance must be handled using a fork lift truck or a pallet truck with suitable forks (fork length at least equal to 2/3 length of unit).

3.2 POSITIONING

Incorrect positioning can cause damage to the appliance and generate hazardous conditions for personnel. The installer must therefore observe the following general regulations:

- fix to the wall the device through the anti-earthquake/tilting chain located in the device backside, using the supplied plastic dowel Ø 8
- make sure you maintain a minimum of 3 cm from the walls
- the room must be well ventilated
- keep well away from sources of heat
- avoid direct sunlight

Specific positioning procedures

- remove packing material (polyethylene, cardboard box, crate, cage)

Polyethylene is potentially dangerous to children

- remove accessories from inside the unit

Wood basement removal: lean the appliance sideways and unscrew the two self-tapping screws (fig.2); drag the appliance from the back side keeping the basement steady till the 4 feet come out from their suitable housings; lean the unit slightly backwards and remove the basement from the front side.

Use gloves when handling wooden packing materials and the wooden base to protect the hands from splinters

- position the appliance with the help of a spirit level. Adjust the levelling feet on the metal base of the unit if necessary (Fig.3)
- remove the protective PVC film from the external surfaces of the unit
- position the shelf (drawer or basket) runners in the holes in the uprights (Fig.4)
- insert the shelves (drawers or baskets) in the runners
- insert the condensate collection tray (in the models where necessary) in the relevant runners located beneath the unit

3.3 WIRING AND ELECTRICAL HOOK-UP

The electrical plant and electrical hook-up operations must be performed by a qualified electrician

For safety reasons adhere to the following indications:

- check that the electrical plant is suitably sized for the absorbed power of the unit
- if the electrical socket and the plug on the appliance power cord are incompatible, change the plug with a suitable component, ensuring the replacement part is of the approved type
- do not use reductions or multi-way adapters (Fig.7)
- check the impedance value of the power supply; the impedance value of the power supply should not exceed 0,075 ohm

It is important to connect the appliance correctly to an efficient earth system executed in compliance with the relevant legislation.

3.4 SET-UP OPERATIONS

To avoid errors and accidents, perform a series of checks for possible damage sustained during transport, installation and hook-up operations before starting up the unit.

Preliminary Checks

- check the condition of the power cord (no cuts or chaffing)
- check that the feet (or rollers), door hinges and shelf supports are stable
- check the condition of internal and external components (pipelines, heat exchanger elements, fans, electrical components, etc.), check also that all parts are firmly fixed into position
- check that the door seals, shelves and drawers are not damaged (broken or scratched) and that the doors close and are sealed properly

The user must also observe the following instructions to obtain the best operation from the appliance:

Indications for Optimal Duty

- do not block the motor compartment air vents
- do not store warm products
- arrange the products on suitable shelves, drawers or baskets. Do not place products directly on the base or against the walls, doors or fixed guards of the unit
- make sure doors are kept closed
- keep the defrost water drain outlet clear
- limit the frequency and duration of opening; each time the door is opened the internal temperature will alter
- load new material at ambient temperature gradually to allow correct refrigeration
- perform routine maintenance regularly (see section 5).

3.5 RE-INSTALLATION

Observe the following procedure:

- switch off the appliance from the main switch
- disconnect the power cord from the electrical outlet
- handle the appliance in accordance with the instructions in heading 3.1
- follow the instructions in headings 3.2 and 3.3 for positioning and hook-ups in the new location

3.6 SCRAPPING AND DISPOSAL

Scrapping and disposal of the appliance must be carried out in full observance of established legislation in your country.

Section 4 OPERATION

4.1 APPLICATIONS AND INTENDED USE

4.1.1 Intended Use and Permitted Use

The appliance is designed and built for preserving and storing medical products on medical/scientific premises.

The appliance is not suitable for the storage of blood or hemoderivates to be reinfused or transfused in the human body.

4.1.2 Improper and Unauthorized Use

1) treatment of products that require constant monitoring and may endure alteration of its characteristics in the case of temperature changes or interruption of refrigeration. For example:

- foodstuff in general
- blood or hemoderivates to be reinfused or transfused in the human body

2) use in places subject to explosive atmosphere

All uses except authorized uses of the appliance shall be construed as "improper use" for which the manufacturer declines all responsibility.

4.2 SAFETY AND ACCIDENT PREVENTION

The appliance embodies various features designed to assure the safety and protect the health of the user. The following list describes the protections adopted against mechanical risks:

- **stability:** the appliance is designed and built so that even with the shelves, drawers or baskets fully extracted in the intended conditions of operation it will remain stable so that it can be used with no risk of tipping, falling or sudden movement
- **surfaces, edges, corners:** accessible parts of the appliance have no sharp corners, sharp edges or rough surfaces that could cause injury
- **moving parts:** moving parts of the unit are designed, built and configured to avoid risk. Moving parts are protected by

fixed guards to prevent accidental contact that could result in injury

Measures adopted for protection against additional risks:

- **electrical power:** the appliance is designed, built and fitted out with the aim of preventing the risk of electric shock in compliance with established safety legislation
- **noise:** the appliance is designed and built to reduce risks related to the emission of airborne noise to a minimum

4.3 SAFETY DATAPLATES AND GUARDS

It is strictly forbidden (Fig.8):

- to tamper with or remove the evaporator cover that protects the user from the risk of cutting on the heat exchanger fins
- to remove the dataplate fixed to the inside edge of the motor housing showing technical specifications (1) and earth connection warning (2)
- to remove the dataplates on the evaporator unit cover near the electrical wiring inside the motor housing which warn the user to disconnect electrical power before working on appliance (3)
- to remove the dataplate fixed inside the motor compartment indicating earthing (4)
- to remove the data tag fixed to the power cord showing the type of power supply (5)

The manufacturer declines all responsibility for safety of the appliance if the above recommendations are not observed.

4.4 OPERATING LIMITS

The appliance is designed and built to work under the following environmental conditions:

SERIES	CLIMATIC CLASS	AMBIENT HUMIDITY
MPR	SN - ST	60 / 70%
MPRR	SN - ST	
LR	T	
LCRR	T	
BLFG	SN - ST	
LFG	SN	
LF	SN	
LDF	SN	
PDF	SN	
LCRF	SN	
CI	SN	
PI	SN	

Internationally, the climatic classes recognized for refrigerators and freezers are:

- **SN or extended temperate class** (ambient temperature from +10°C to +32°C);
- **N or temperate class** (ambient temperature from +16°C to +32°C);
- **ST or subtropical class** (ambient temperature from +18°C to +38°C);
- **T or tropical class** (ambient temperature from +18°C to +43°C).

The standard power supply must be 230V +/- 10% 50Hz.

Section 5 ROUTINE AND PROGRAMMED MAINTENANCE

The information in this section regards the user, or other non-specialized personnel, and the routine maintenance technician.

5.1 BASIC SAFETY REGULATIONS

We summarize the safety regulations already shown in heading 1.5 to ensure that the user or maintenance technician can perform the work in conditions of total safety:

- do not touch the unit with wet hands and/or feet
- do not use the appliance with bare feet
- do not insert screwdrivers or other pointed objects between guards or moving parts of the appliance

- do not pull the power cord to disconnect the appliance from the electrical mains
- before performing any cleaning or maintenance on the appliance disconnect it from the electrical mains by switching of the main switch and extracting the plug

5.1.1 Prohibited: Removal of Guards and Safety Devices

It is strictly forbidden to remove guards or safety devices when performing routine maintenance work. The manufacturer disclaims all liability that may arise if this regulation is not observed.

5.1.2 Indications on Emergency Measures in Case of Fire

- disconnect the appliance from the electrical power socket or switch off the master switch on the electrical mains line
- do not use water to douse fires
- use powder or foam extinguishers

5.2 CLEANING THE APPLIANCE

In view of the usage destination of the appliance, it is important to keep it clean for reasons of hygiene and health. The appliance is thoroughly cleaned in our factory before delivery. We recommend, however, to clean the interior of the appliance before use. Before cleaning the appliance make sure the power cord is disconnected.

5.2.1 Cleaning the Interior and Exterior of the Appliance

- cleaning products: water and non-abrasive neutral detergent. **DO NOT USE SOLVENT OR THINNERS**
- cleaning method: use a cloth or sponge soaked in a suitable cleaning product to clean the interior and exterior parts of the cabinet
- sanitation: do not use substances that could alter the stored products
- rinsing: use a cloth or sponge soaked in clean water. **DO NOT USE WATER JETS**
- frequency: once a week or at different intervals in accordance with the type of product conserved.

5.2.2 Cleaning the Condenser

The condenser will work less efficiently if it is obstructed with foreign material so it must be cleaned once a month. Before cleaning the condenser switch off the appliance, disconnect the power cord and proceed as follows:

Motor at base - open the control panel by slackening the screw positioned in the top part and pull it up

Motor at top - use a step ladder for direct access to the condenser located at the top of the appliance.

Use an air jet or dry brush and, working with up and down movements (Fig.9), remove any dust or fluff that has deposited on the heat exchanger fins. If there are greasy deposits on the fins use a brush soaked in benzene or alcohol. After cleaning, start the appliance (with bottom motor models, remember to close control panel).

During this operation use the following personal safety measures: safety glasses, respirator mask, chemical resistant gloves (benzene - alcohol).

5.3 PERIODIC CHECKS

The following areas of the appliance or component assemblies require periodic checking (anniversaries):

- condition and efficiency of the door sealing elements
- condition of shelves in contact with the products
- condition of hinges and correct fixing of the doors
- condition of electrical cables and electrical parts

5.4 PRECAUTIONARY MEASURES FOR PROLONGED DISUSE

If the appliance is to remain unused for more than 15 days proceed as follows:

- switch off the appliance and disconnect it from the electrical supply
- clean the interior of the cabinet, shelves, drawers, baskets, runners and supports, paying special attention to critical areas such as articulations and magnetic sealing strips in accordance with the indications in heading 5.2.
- leave doors slightly open to prevent accumulation of residual humidity

5.5 PREVENTIVE MAINTENANCE

5.5.1 Start-up after Prolonged Disuse

Before starting the appliance after prolonged disuse perform preventive maintenance. Clean the unit thoroughly as described in heading 5.2.

5.5.2 Checking Warning and Control Devices

Check the correct running of the controls according to what is reported in the "Instruction and Maintenance Manual" enclosed. We recommend you to take out a service or maintenance contract with your dealer covering:

- cleaning of the condenser
- keeping a check on the refrigerant charge
- checking complete cycle operation
- electrical safety

Section 6 SPECIAL MAINTENANCE AND REPAIRS

All maintenance work not described in the previous sections must be considered "Special Maintenance".

Special maintenance interventions and repairs are to be performed exclusively by specialized technicians authorized by the manufacturer.

The manufacturer declines all liability in the case of work performed by the user or unauthorized persons, or if non-original spare parts are fitted to the appliance.

Section 7 DIAGNOSTICS

The following problems may occur as indicated in the table below:

FAULT	POSSIBLE CAUSE	ACTION
Appliance does not switch on	switch set to - O - (or OFF) power failure other	set switch to - I - (press the button ON/OFF) check plug, socket, fuses, electrical line contact service department
Refrigeration unit does not start	set temperature has been reached defrosting cycle in progress control panel fault other	set new temperature wait for cycle end, switch off and then on again contact service department contact service department
Refrigeration unit runs constantly but does not reach set temperature	room temperature too high condenser dirty refrigerant needs recharging condenser fan not running door seals are not efficient evaporator is coated with ice	provide better ventilation of room clean condenser contact service department contact service department check seals / make sure appliance contents are not obstructing doors defrost manually
Refrigeration unit does not cut out at set temperature	control panel fault temperature sensor defective	contact service department contact service department
Build up of ice on evaporator	improper use control panel fault defrost sensor defective	see headings 3.4 contact service department contact service department
Water or ice deposits in the drip tray	drain outlet is clogged appliance not level	clean drain and drain outlet see heading 3.2

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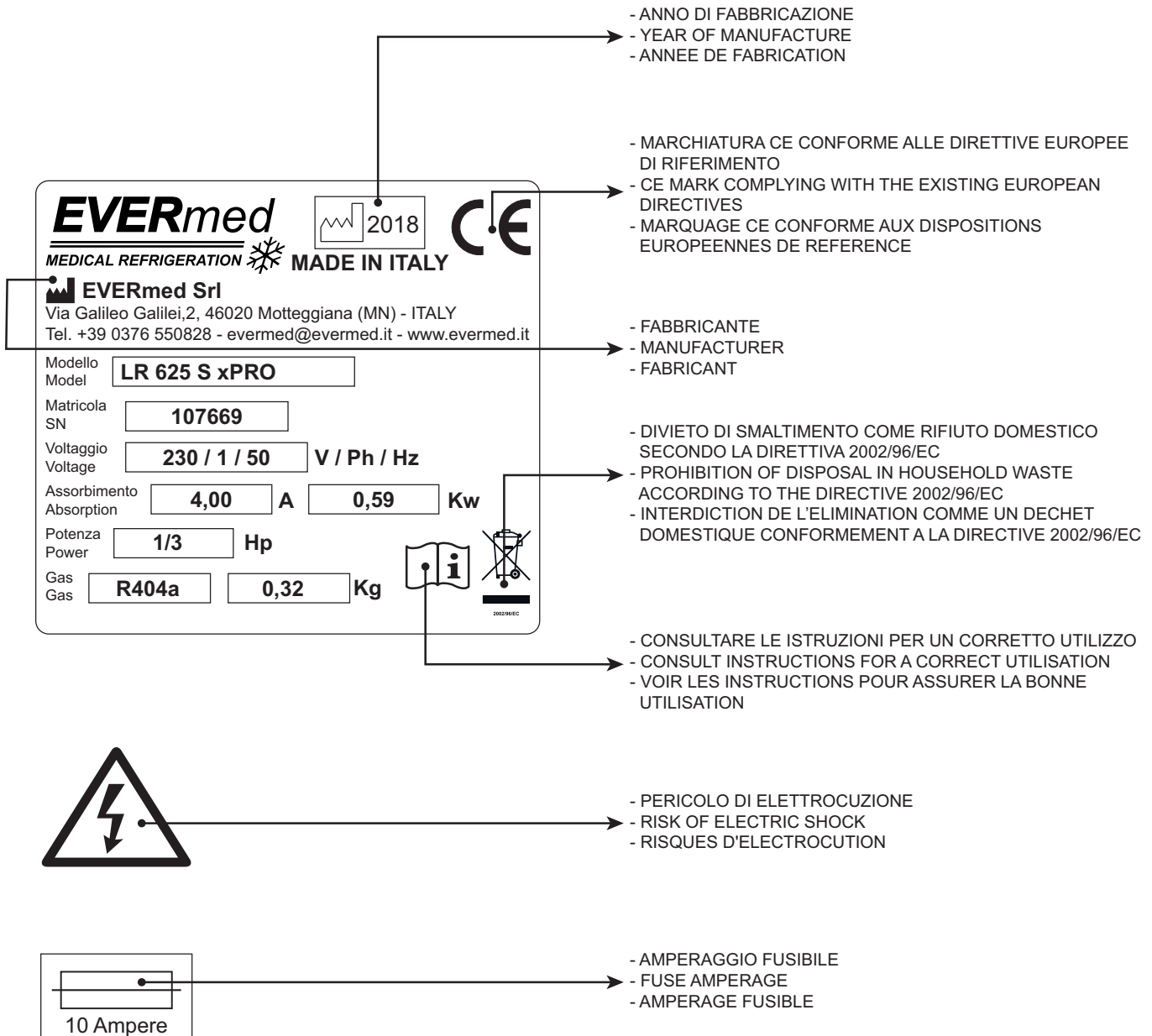
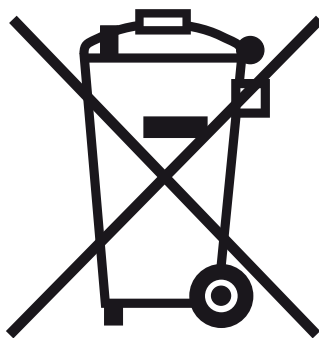
SYMBOLS:

Over the appliance you may find the following labels with the icons below described.

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ENGLISH

The symbol  means that this product should not be treated as a household waste.

To prevent potential negative consequences for the environment and health, be sure this product is correctly disposed of and recycled.

Further information on the disposal and recycling of this product, please contact your Distributor, after Sales Service or the waste Treatment Service.

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EVERmed Srl: Via Galileo Galilei, n° 2 - 46020 Motteggiana (MN) ITALY